# Safety Related Information

for the Thermo Scientific Medifuge Centrifuge

## Intended Use

This centrifuge is intended to be used as a laboratory equipment to separate sample mixtures of different densities.

This centrifuge can be used as an In-Vitro diagnostics device (Directive 98/79/EC), if used together with IVD tubes, to separate blood into its components such as serum and plasma for further clinical diagnostic analysis.

The centrifuge has to be operated by a trained individual such as a clinical laboratory technologist or a trained laboratory technician.

## Precautions

### Signal Words and Colors

#### WARNING

Indicates a hazardous situation that, if not avoided, could result in death or serious injury.

#### CAUTION

Indicates a hazardous situation that, if not avoided, could result in minor or moderate injury (e.g. sample loss).

## Set Up Conditions

As a safety zone maintain a clear radius of at least 30 cm around the centrifuge. Do not place any dangerous substances within this security zone.

### CAUTION

- Implement measures which ensure that no one can approach the centrifuge for longer than absolutely necessary while it is running.
- Do not use a damaged rotor. Replace the rotor, if it was dropped.
- Not following the set up conditions can result in minor or serious injuries.

### WARNING

It is the obligation of the operator to make sure, that protective clothing is used. Mind the "Laboratory Biosafety Manual" of the World Health Organization (WHO) and the regulations in your country.

## Hazardous Substances

### WARNING

- Especially when working with corrosive samples (salt solutions, acids, bases), the accessory parts and vessel have to be cleaned thoroughly.
- Do not centrifuge explosive or flammable materials or substances.
- The centrifuge is not inert or protected against explosion. Never use the centrifuge in an explosion-prone environment.
- Do not centrifuge toxic or radioactive materials or any pathogenic micro-organisms without suitable safety precautions.

## Operation

### WARNING

- Never use the centrifuge if parts of its cover panels are damaged or missing.

- Do not move the centrifuge while it is running.
- Do not place anything on top of the centrifuge during a run.

### CAUTION

- Do not lean on the centrifuge.
- Never open the centrifuge door until the rotor has come to a complete stop and this has been confirmed in the display.
- The emergency door release may be used in emergencies only to recover the samples from the centrifuge, e.g. during a power failure.

## Explanation of Symbols

This symbol refers to general hazards.

This symbol refers to biological hazards.

Observe the information contained in the instruction manual to keep yourself and your environment safe.

This symbol refers to physical hazards from hot surfaces.

This symbol refers to information on hazards, described within the manual.

This symbol demands to disconnect mains before transporting or servicing the centrifuge.

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