


**Thermo Scientific B3 Mini Gel System  
Casting with Owl's Gel Casting System**


**1. Place UVT gel tray in buffer chamber in the casting position. Gaskets will form a seal against the walls of the chamber.**



**2. Pour warm (<math><60^\circ</math>) agarose onto tray and set combs in the desired comb slot(s).**

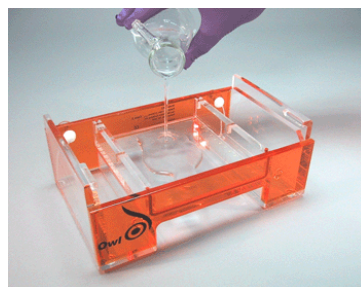

**3. Once solidified, turn the tray 90 degrees to the running position, remove combs, add buffer, load samples and run the gel.**

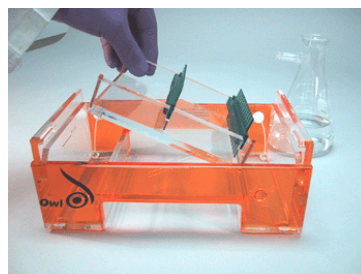

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

>> For more information, visit: [www.thermoscientific.com/owlsci](http://www.thermoscientific.com/owlsci)