



**Thermo Scientific Owl D4 System
Gel Casting Using the Owl Model D4**

1. Place the UVT gel tray(s) inn the multiple external gel caster. Level the gel caster. The gasketed sides of UVT gel tray form a leak proof seal along the side of the gel tray.

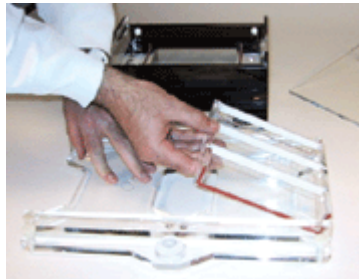

2. Prepare and pour agarose solution. Allow agarose to cool then place comb(s) into slot(s) provided on walls of UVT gel tray.

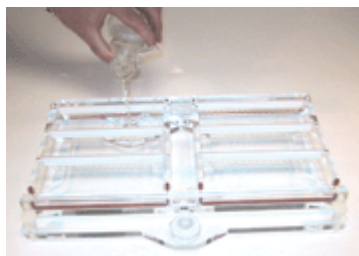

3. Once the gel has solidified, remove the comb(s), place the UVT gel tray(s) into the buffer chamber, load samples, add buffer and run the gel.



© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

>> For more information, visit: www.thermoscientific.com/owlsci