


**Thermo Scientific Owl JGC-4 Gel Casting System**  
**Gel Casting Using the Owl Model JGC-4**


**1. Assemble the gel cassette (s) and place into a pouch.**

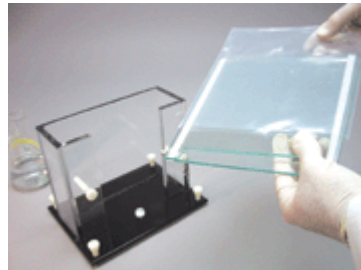

**2. Tighten the cassette in the caster.**



**3. Pour the acrylamide solution onto the pouch. Once the gel is polymerized, remove the cassette and place in a running device.**

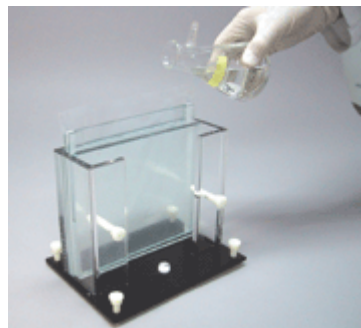

© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

>> For more information, visit: [www.thermoscientific.com/owlsci](http://www.thermoscientific.com/owlsci)