

**Using the VEP-3 Electroblotting System**

- 1. Stack gel and membrane between foam pads and blotting filter paper, place into the cassette and place clip onto the cassette**
- 2. Slide the cassette into the tank**
- 3. The tank is designed so there is no chance of setting the cassette(s) in backwards**
- 4. If temperature regulation is required, attach tubing to the hose connectors on the back of the tank**
- 5. The electroblotting system may be hooked up to tap water or a recirculator for cooling**

*© 2012 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.*

>> For more information, visit: [www.thermoscientific.com/owlsci](http://www.thermoscientific.com/owlsci)